

## Fyrite Reading Instructions:

**Note: Be sure to hold canister away from face. Do not press button with canister inverted. To ensure accuracy, it is important to always follow the same procedure when taking a Fyrite reading.**

~ Press button on top of Fyrite canister to release CO<sub>2</sub> concentration. Tip canister to the side to ensure all fluid is released from top of canister.

~ Loosen screw on slide scale & align top of fluid with zero on the scale. Tighten screw.

~ Connect hose & bulb to unit being tested.

~ Place hose directly over button on top of canister & press firmly.

~ With button depressed, squeeze bulb 27 times. On the last squeeze, with bulb still deflated, release hose from button.

~ Turn Fyrite canister upside down 3 times, each time allowing all fluid to flow to the opposite end of the canister.

~ Tip canister slightly to ensure all fluid has been released from top of canister.

~ Read CO<sub>2</sub> concentration in %.

**Note: Your Fyrite indicator will come with a complete set of detailed instructions which should be followed carefully. The fluid used inside this Fyrite instrument is poisonous & corrosive & must not be taken internally. In event of a spill or accidental body contact with the Fyrite fluid, follow instructions on the refill bottle carefully.**